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The **VACUUM IMPREGNATION UNIT** is a small, practical device for embedding or impregnating porous metallographic samples in a vacuum. The device, which was developed by SCAN-DIA, is easy to use and facilitates perfect embedding.

The VACUUM IMPREGNATION UNIT consists of a plastic housing with built-in low vibration electric vacuum pump. On the housing stands a glass vacuum chamber (inside diameter 100 mm x 60 mm high), which offers an optimal view of the embedding process. The device is also equipped with a vacuum gauge and regulator valve. The level of the vacuum can therefore be individually adjusted and read off immediately. There are also a cover plate, a rubber seal, silicon-rubber funnels, glass stoppers and glass inlet-tubes.

The process is very simple. First place the sample in a SCANDIFORM and then put them both into the glass vacuum chamber. Close the chamber with the cover plate which has the funnel together with the glass stopper on top. Switch the device on and wait until the desired vacuum is obtained. Under vacuum pour the mixed coldmounting compound on to the sample via the funnel. The ventilation which follows allows the mounting compound to penetrate the pores and cracks in the sample. Please note that only pores and cracks which are connected to the outer surface are filled.

All cold mounting compounds can be used. Best results are achieved with products which do not out-gas in a vacuum of approximately 0,6 bar, have low viscosity and a long curing time. The cold mounting compounds SCANDIPILEX and SCANDIPLAST are especially recommended.

A larger glass vacuum container is available as an option.

Machine:

Item-no. 9117 VACUUM IMPREGNATION UNIT including 3 Silicon rubber funnels, 10 Glass stoppers, 100 Glass inlet tubes and 1 Cover plate.

parts:

Item-no. 91171 Glass stoppers, 25 pieces
Item-no. 91172 Glass inlet tube, 100 pieces
Item-no. 91173 Silicon rubber funnel, 1 piece
Item-no. 91174 Cover plate, 1 piece
Item-no. 91175 Glass vacuum chamber, 100 mm dia. x 60 mm height, 1 piece
Item-no. 91176 Rubber gasket, 1 piece

option:

Item-no. 91177 Glass vacuum chamber, 140 mm dia. x 60 mm height, including larger cover plate and rubber gasket

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